

PuroMill™ Pharmaceutical Milling Media

Solubility enhancement through advanced
polymeric milling media technology



PuroLite®
An Ecolab Company



Puro-lite®

An Ecolab Company

About Puro-lite

Puro-lite is a leading manufacturer of ion exchange, catalyst, adsorbent and specialty resins. With global headquarters in the United States of America, Puro-lite is the only company that focuses 100% of its resources on the development and production of resin technology.

Responding to our customers' needs, Puro-lite has the widest variety of products and the industry's largest technical sales force. Globally, we have strategically located research and development centers and application laboratories. Our ISO 9001 certified manufacturing facilities in the USA, United Kingdom, Romania and China combined with more than 40 sales offices in 30 countries ensure complete worldwide coverage.

Puro-lite has been part of Ecolab since 2021. A trusted partner at nearly three million commercial customer locations, Ecolab (ECL) is the global leader in water, hygiene and infection prevention solutions and services. Ecolab delivers comprehensive solutions, data-driven insights and personalized service to advance food safety, maintain clean and safe environments, optimize water and energy use, and improve operational efficiencies and sustainability for customers in the food, healthcare, hospitality and industrial markets in more than 170 countries around the world.



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
PREMIER PRODUCTS

The quality and consistency of our products are fundamental to our performance. Throughout all Puro-lite plants, production is carefully controlled to ensure that our products meet the most stringent criteria, regardless of where they are produced.



RELIABLE SERVICE

We are technical experts and problem solvers. Reliable and well-trained, we understand the urgency required to keep businesses operating smoothly. Puro-lite employs the largest technical sales team in the industry.



INNOVATIVE SOLUTIONS

Our continued investment in research and development means we are always perfecting and discovering innovative uses for ion exchange resins and adsorbents. We strive to make the impossible possible.

PuroMill Pharmaceutical Milling Media

Contents

What is PuroMill Pharmaceutical?	3
Universal Formulation Approach with Pure Drug Nanoparticles	4
Typical Chemical and Physical Characteristics	5
Mill with Confidence	8

What is PuroMill Pharmaceutical?

PuroMill pharmaceutical-grade milling media is an advanced polymer-based bead-form milling media that is extremely durable, reusable and chemically inert. PuroMill enables the preparation of high-purity nanoparticles for the development of small molecule pharmaceutical formulations using conventional nanoparticle milling technologies. PuroMill pharmaceutical-grade milling media is manufactured in conformance with cGMP guidelines under stringent specifications using advanced copolymer synthesis and purification technologies that assure the highest attainable drug product quality.

PuroMill is an essential element of nanoparticle milling processes used to improve the solubility and bioavailability of poorly soluble small molecules. PuroMill media is ideal for pharmaceutical companies, CMOs and CDMOs engaging in API and formulation development of oral, injectable and inhalation dosage forms based on liquid or solid formulations. Additionally, PuroMill also effectively enables efficient isolation of macromolecules through biological cell disruption.

Improve Bioavailability of Drug Compounds Without Process-Related Contamination

- Highly purified polymeric milling media
- Produced in a cGMP certified, FDA inspected facility
- Ideal for pharmaceutical/biopharmaceutical drug formulations
- Creates drug nanoparticles < 100 nm to enhance absorption of drugs with poor solubility
- Enables efficient biological cell disruption without milling-related contamination
- Smooth, non-porous, non-adsorptive micro surface minimizes contamination
- Exceptional wear resistance minimizes particulate contamination from media attrition and equipment abrasion
- Non-reactive and biologically inert
- Autoclavable/steam sterilizable
- Low density; reduces frictional heat to prevent degradation of heat-sensitive APIs
- Available in sizes from 50 µm–1,000 µm for optimized milling performance of drug formulations

Universal Formulation Approach with Pure Drug Nanoparticles

Because of the high percentage of poorly soluble compounds in the development pipeline, formulation teams have significant challenges advancing drug products for clinical evaluation. PuroMill pharmaceutical-grade, high-performance milling media reduces API particles to the nano scale.

Pure drug nanoparticles have increased surface area, faster dissolution and improved solubility, which can enhance bioavailability and enable customization of pharmacokinetic profiles. The results include improved drug efficacy, safety and patient compliance.

PuroMill pharmaceutical advanced high-energy media technology minimizes the risk of contamination due to leaching as well as from attrition of milling media and milling equipment surfaces. With exceptional wear resistance, unmatched purity and milling efficiency, nanoparticles of API compounds can be successfully created with previously unattainable levels of product quality and purity.

PuroMill enables production of pure drug nanoparticles, stabilized by low concentrations of GRAS stabilizers to enable high drug loaded formulations. Pure drug nanoparticle formulations can accelerate in-vivo drug dissolution and improve saturation solubility to maximize the rate and extent of absorption. PuroMill can also help achieve ultra-fine nanoparticle formulations suitable for IV, IM and SC injections.

Typical Chemical and Physical Characteristics

PuroMill Pharmaceutical PM3000

Characteristic	Test Method	Result
< 150 microns	PITM002E	0.01% (max.)
250–350 microns	PITM002D	95% (min.)
Volume Median Diameter	PITM002E	250–350 µm
Residual Monomers	PITM086B	
Styrene		50 ppm (max.)
Divinylbenzene		50 ppm (max.)
Ethylvinylbenzene		50 ppm (max.)
Residual Solvent - Methanol	PITM057F	100 ppm (max.)
Heavy Metals	PITM017C	10 ppm (max.)
Microbial Limits	PITM055A	
Total Aerobes		< 10 CFU/g (max.)
Total Yeasts		< 10 CFU/g (max.)
Total Molds		< 10 CFU/g (max.)
Escherichia Coli		Negative/10g
Salmonella Species		Negative/10g
Pseudomonas Aeruginosa		Negative/10g
Staphylococcus Aureus		Negative/10g
Bacterial Endotoxins	PITM055B	0.5 EU/g (max.)
Appearance	PITM001	Non-aggregated spherical beads, white to off-white in color

PuroMill Pharmaceutical PM5000


Characteristic	Test Method	Result
< 200 microns	PITM002E	0.02% (max.)
425–575 microns	PITM002D	95% (min.)
Volume Median Diameter	PITM002E	450–550 µm
Residual Monomers	PITM086B	
Styrene		50 ppm (max.)
Divinylbenzene		50 ppm (max.)
Ethylvinylbenzene		50 ppm (max.)
Residual Solvent - Methanol	PITM057F	100 ppm (max.)
Heavy Metals	PITM017C	10 ppm (max.)
Microbial Limits	PITM055A	
Total Aerobes		< 10 CFU/g (max.)
Total Yeasts		< 10 CFU/g (max.)
Total Molds		< 10 CFU/g (max.)
Escherichia Coli		Negative/10g
Salmonella Species		Negative/10g
Pseudomonas Aeruginosa		Negative/10g
Staphylococcus Aureus		Negative/10g
Bacterial Endotoxins	PITM055B	0.5 EU/g (max.)
Appearance	PITM001	Non-aggregated spherical beads, white to off-white in color

PuroMill Pharmaceutical PM7000

Characteristic	Test Method	Result
< 250 microns	PITM002E	0.02% (max.)
600–800 microns	PITM002D	95% (min.)
Volume Median Diameter	PITM002E	650–750 µm
Residual Monomers	PITM086B	
Styrene		50 ppm (max.)
Divinylbenzene		50 ppm (max.)
Ethylvinylbenzene		50 ppm (max.)
Residual Solvent - Methanol	PITM057F	100 ppm (max.)
Heavy Metals	PITM017C	10 ppm (max.)
Microbial Limits	PITM055A	
Total Aerobes		< 10 CFU/g (max.)
Total Yeasts		< 10 CFU/g (max.)
Total Molds		< 10 CFU/g (max.)
Escherichia Coli		Negative/10g
Salmonella Species		Negative/10g
Pseudomonas Aeruginosa		Negative/10g
Staphylococcus Aureus		Negative/10g
Bacterial Endotoxins	PITM055B	0.5 EU/g (max.)
Appearance	PITM001	Non-aggregated spherical beads, white to off-white in color

FIGURE 1

The smooth, non-porous, non-adsorptive micro surface of PuroMill minimizes contamination while the highly consistent, monodisperse particle size prevents screen binding and ensures reproducible milling performance.

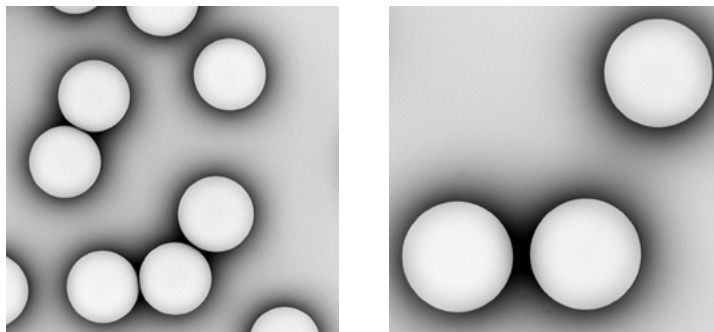

Mill with Confidence

PuroMill is a pharmaceutical-grade, cGMP compliant contact material.

PuroMill pharmaceutical styrene-divinylbenzene copolymer based milling media is for high-energy wet media milling in pharmaceutical API and formulation development applications and meets exacting standards for new generations of nanomaterials processing.

Highly purified to eliminate leaching and extraction of soluble materials (including solvents and monomers), PuroMill is non-reactive and biologically inert to prevent undesirable chemical reactions and composition changes. The polymer beads generate less heat than traditional media to ensure the processed molecule remains thermodynamically stable and structurally unaltered. Additionally, the low media density (1.07 g/cc) enables high media loads and agitation speeds to maximize particle size reduction without contamination – while reducing scale-up inefficiencies associated with media centrifugation of more dense media. Plus, the monodisperse (uniform) particle size distribution and low fines composition prevents screen binding and enables easy separation from the milled compound.

In addition to use in conventional media mills, PuroMill is compatible with novel nanoparticle milling and cell disruption processes such as high-speed dispersers, homogenizers and rotor-stators.

- The only commercially available cGMP milling media for pharmaceutical and biotechnology applications
- Can be used in AISI 316 stainless steel or other suitable metal alloy mills for sanitary and aseptic applications, eliminating the complexity and cost associated with ceramic media mill equipment designs
- Conforms to stringent USP <61> and USP <85> specifications for microbial limits and endotoxins
- Complies with strict USP <233> specifications for elemental impurities
- Drug Master File registered with the U.S. FDA



Algeria
Australia
Bahrain
Brazil
Canada
China
Czech Republic
France
Germany

India
Indonesia
Israel
Italy
Japan
Jordan
Kazakhstan
Korea
Malaysia

Mexico
Morocco
New Zealand
Poland
Romania
Russia
Singapore
Slovak Republic
South Africa

Spain
Taiwan
Tunisia
Türkiye
UK
Ukraine
USA
Uzbekistan



Americas

Purolite
2201 Renaissance Blvd.
King of Prussia, PA 19406
T +1 800 343 1500
T +1 610 668 9090
F +1 800 260 1065
americas@purolite.com

EMEA

Purolite Ltd.
Unit D
Llantrisant Business Park
Llantrisant, Wales, UK
CF72 8LF
T +44 1443 229334
F +44 1443 227073
emea@purolite.com

FSU

Purolite Ltd.
Office 6-1
36 Lyusinovskaya Str.
Moscow, Russia
115093
T +7 495 363 5056
F +7 495 564 8121
fsu@purolite.com

Asia Pacific

Purolite China Co. Ltd.
Room 707, C Section
Huanglong Century Plaza
No.3 Hangda Road
Hangzhou, Zhejiang, China 310007
T +86 571 876 31382
F +86 571 876 31385
asiapacific@purolite.com

Purolite, the leading manufacturer of quality ion exchange, catalyst, adsorbent and specialty high-performance resins, is the only company that focuses 100% of its resources on the development and production of resin technology.

We're ready to solve your process challenges. For further information on Purolite products and services, visit www.purolite.com or contact your nearest Technical Sales Office.



www.purolite.com

©2022 Purolite
All rights reserved.
P-000130-100PP-1022-R5-PCO

The statements, technical information and recommendations contained herein are believed to be accurate as of the date hereof. Since the conditions and methods of use of the product and of the information referred to herein are beyond our control, Purolite expressly disclaims any and all liability as to any results obtained or arising from any use of the product or reliance on such information; NO WARRANTY OF FITNESS FOR ANY PARTICULAR PURPOSE, WARRANTY OF MERCHANTABILITY OR ANY OTHER WARRANTY, EXPRESSED OR IMPLIED, IS MADE CONCERNING THE GOODS DESCRIBED OR THE INFORMATION PROVIDED HEREIN. The information provided herein relates only to the specific product designated and may not be applicable when such product is used in combination with other materials or in any process. Nothing contained herein constitutes a license to practice under any patent and it should not be construed as an inducement to infringe any patent and the user is advised to take appropriate steps to be sure that any proposed use of the product will not result in patent infringement.